
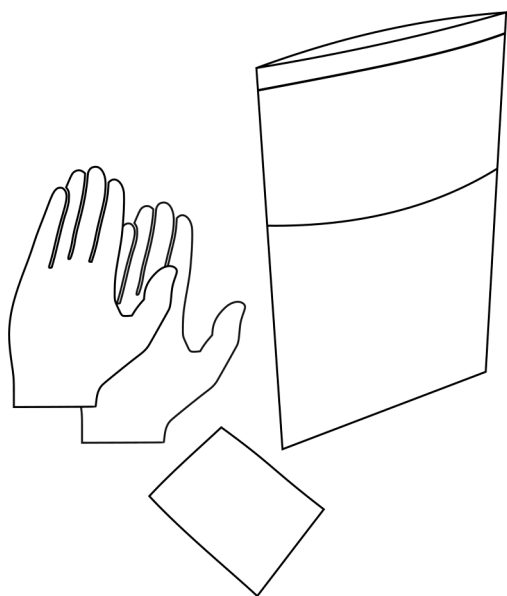


# Hy Con

MECCONTI

## ENVIRONMENTAL SURFACE SAMPLING

For Samples  
Collection from  
Not Disinfected  
Surfaces and  
Carcasses of  
Slaughter Animals  
(according to ISO  
17604: 2015)

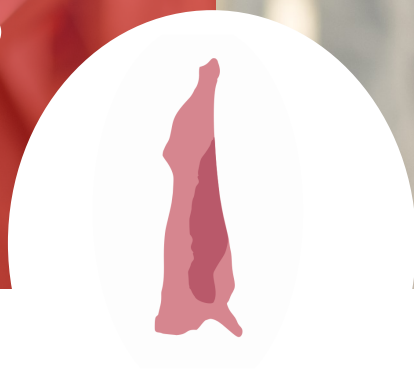

## HY CON CHECK S/BPW PRE-MOISTENED SAMPLING SPONGES SK004

EACH STERILE SET CONTAINS:

SAMPLING MATERIAL - **CELLULOSE SPONGE PRE-MOISTENED WITH BUFFERED PEPTONE WATER**, A PAIR OF **DISPOSABLE GLOVES** & A **SAMPLE BAG WITH A POCKET** FOR THE SAMPLE DESCRIPTION

# HY CON CHECK S/BPW



Commission Regulation (EU) No 2073/2005 of 15 November 2005 impose an obligation on the slaughterhouses to collect samples from carcasses of slaughter animals in accordance with the techniques described in ISO 17604: 2015. The non-destructive method described in the standard is to collect sample from the appropriate carcass locations using an abrasive sponge. Collected material should be returned to the laboratory within 24 hours of receipt (optimally within an hour). After collection and during transportation samples should be stored refrigerated.

## IDEAL FOR

**NON-DISINFECTED SURFACES**

**CARCASSES OF SLAUGHTER ANIMALS**

**QUALITATIVE RESEARCH**

e.g. : the detection of Salmonella spp., or Listeria monocytogenes

- **Convenient and Easy To Use**
- **Ready to Use** - do not require additional moistening before sampling
- **Smart 'Kangaroo' bag** - additional pocket keep the gloves before sampling and can keep a description after
- **Reliable results** - components of the kit do not inhibit the microbial growth, sterile gloves reduce the cross-contamination risk

## *Why Mecconti?*

*„We are an experienced producer of high quality innovative products. Our competitive solutions are always tested in close cooperation with end-users.*



*We are flexible and open to all customer needs.*

*The environment is key for us. We do our best to minimize environmental impact by using biodegradable materials and minimizing waste”.*


### References:

1. ISO 18593:2005 Microbiology of food and animal feeding stuffs - Horizontal methods for sampling techniques from surfaces using contact plates and swabs.
2. ISO 17604:2015-10 Microbiology of the food chain – Carcass sampling for microbiological analysis.
3. Commission Regulation (EC) No 2073/2005 of 15 November 2005 on microbiological criteria for foodstuffs.


# SAMPLE COLLECTION

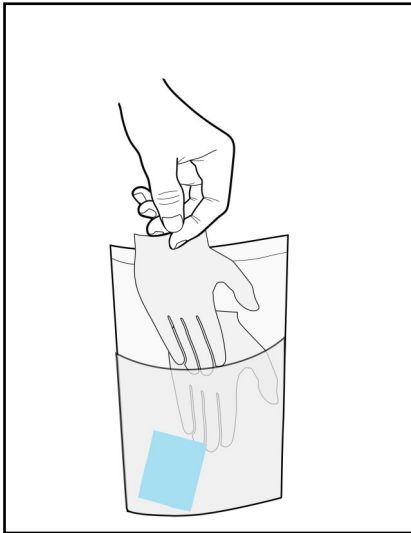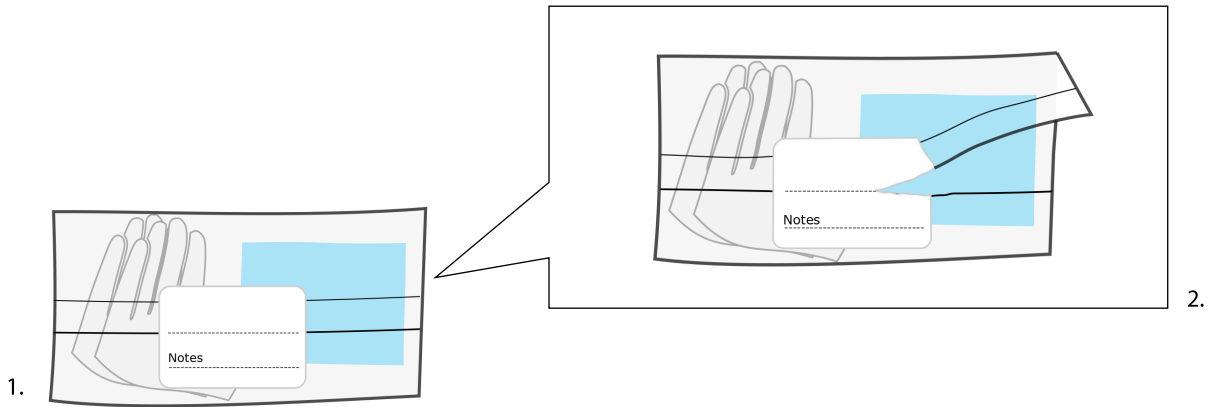


3. remove gloves

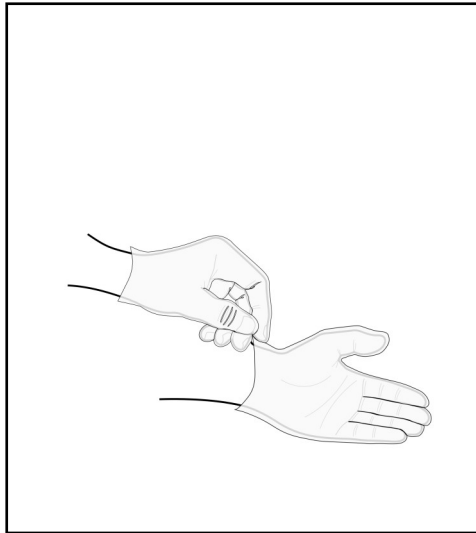


4. put the gloves on

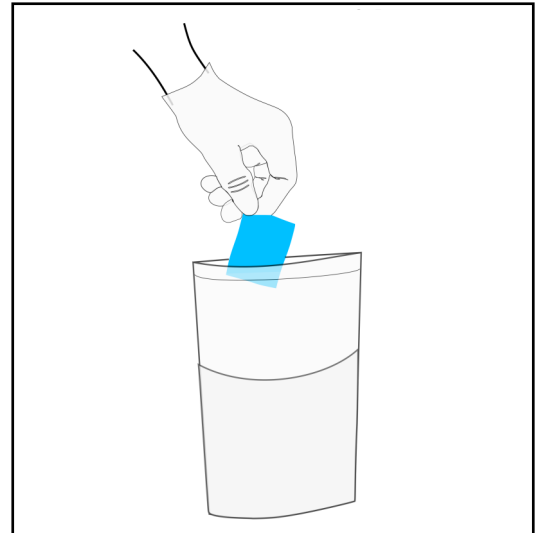


5. take out the sponge

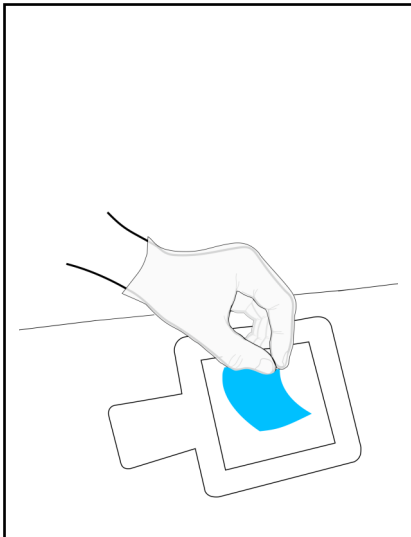


6. swab the defined surface

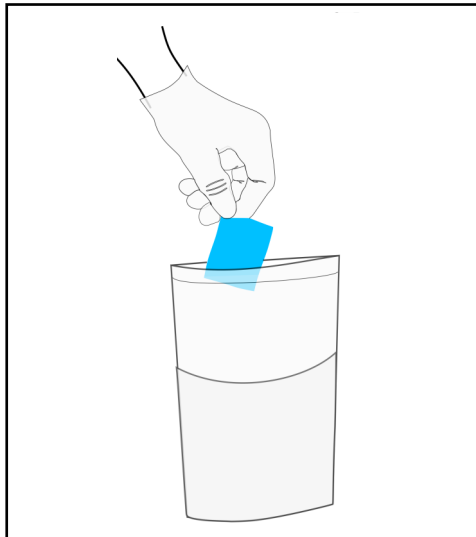

7. return used sponge into sample bag

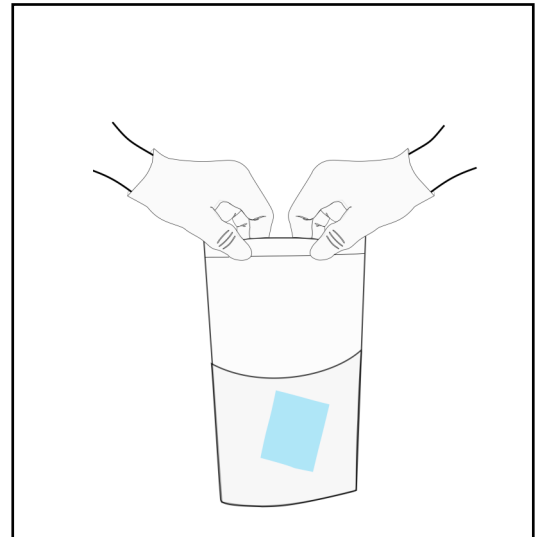

8. close the zip

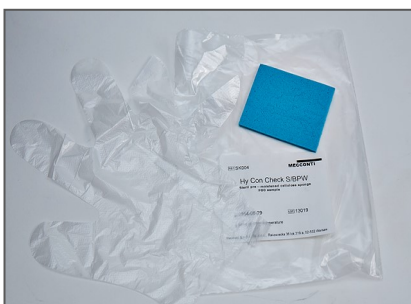

HY CON CHECK S/BPW			
CODE	PRODUCT DESCRIPTION	ORDERING/PACKING	
SK004	each sterile set contains: pre-moistened with 10 ml BPW cellulose sponge to collect environmental samples, 2 gloves and the sample bag with a pocket for the sample description	SK004.0010	1 pack-10 pcs (2 x 5pcs)
		SK004.0250	1 pack-250 pcs (50 x 5pcs)

**MECCONTI S.A.R.L. SP. Z O.O.**

T: +48 22 3974533, F: +48 22 5047213  
sales@mecconti.com, support@mecconti.com  
www.mecconti.com,  
www.facebook.com/Mecconti  
www.linkedin.com/company/mecconti