

Hy Con

MECCONTI

ENVIRONMENTAL SURFACE SAMPLING

A Comprehensive
Solution for
Swabbing from
Wet Surfaces
and Hands.



HY CON CHECK W

DRY SAMPLING WIPES

SK001

EACH **STERILE SET** CONTAINS:

SAMPLING MATERIAL - **VISCOSE CLOTH**

A PAIR OF **DISPOSABLE GLOVES**

SAMPLE BAG WITH A POCKET FOR THE SAMPLE DESCRIPTION





Commission Regulation (EU) No 2073/2005 of 15 November 2005 impose an obligation on the slaughterhouses to collect samples from carcasses of slaughter animals in accordance with the techniques described in ISO 17604: 2015 and to the production facilities suggest the collection of environmental samples in accordance with ISO18593. Samples can be collected using sterile abrasive sponges, wipes and swabs.

VERSATILE APPLICATION e.g. WET SURFACES AND HANDS

RECOMMENDED FOR QUANTITATIVE RESEARCH

IDEAL FOR LARGER SURFACES

- **Convenient and easy to use**
- **Reduced waste**
- **Flexible material** allows swabbing from uneven, complex surfaces
- Possibility to moisten with **preferred buffer** just prior to sample collection
- **Higher recovery** in comparison with the standard used cellulose sponge or cotton swab

Why Mecconti?

„We are an experienced producer of high quality innovative products. Our competitive solutions are always tested in close cooperation with end-users.


We are flexible and open to all customer needs.

The environment is key for us. We do our best to minimize environmental impact by using biodegradable materials and minimizing waste”.


References:

1. ISO 18593:2005 Microbiology of food and animal feeding stuffs - Horizontal methods for sampling techniques from surfaces using contact plates and swabs.
2. ISO 17604:2015-10 Microbiology of the food chain – Carcass sampling for microbiological analysis.
3. Commission Regulation (EC) No 2073/2005 of 15 November 2005 on microbiological criteria for foodstuffs.


SAMPLE COLLECTION

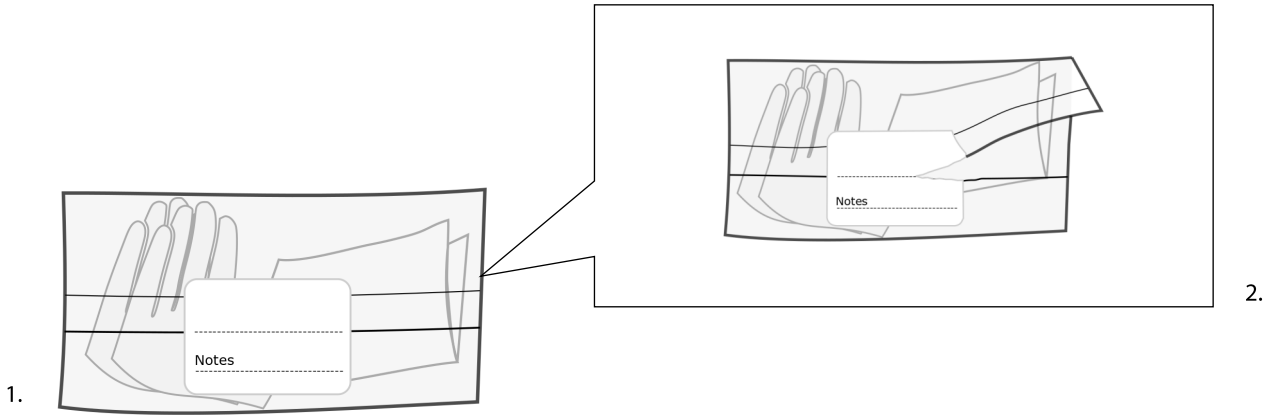


3. remove gloves

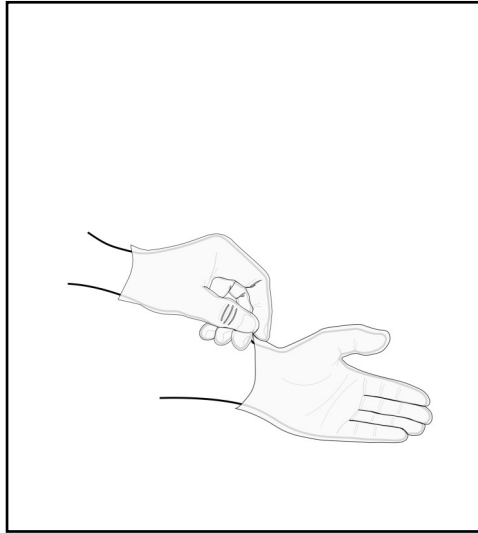


4. put the gloves on

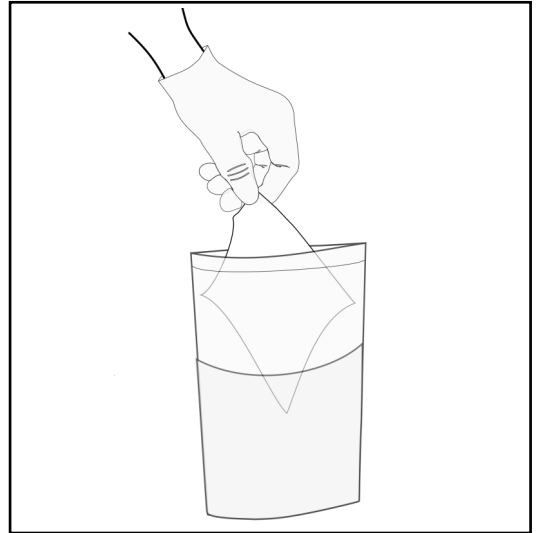


5. take out the wipe

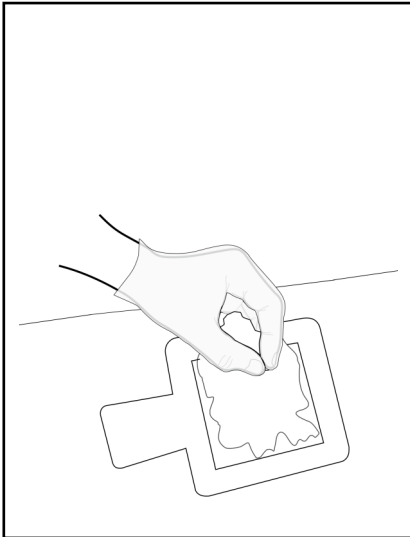

6. wipe the defined surface

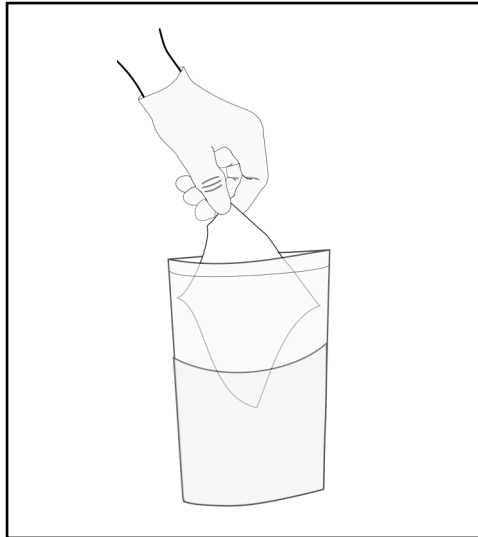

7. return used wipe into sample bag

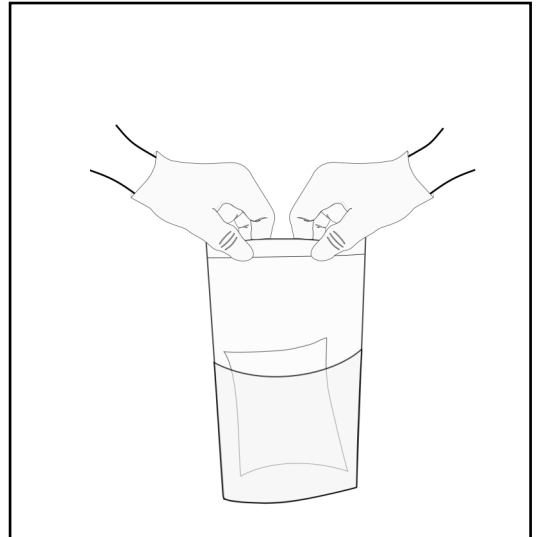

8. close the zip



HY CON CHECK W

CODE	PRODUCT DESCRIPTION	ORDERING/PACKING	
SK001	each sterile set contains: viscose cloth 25x40cm to collect environmental samples, 2 gloves and the sample bag with a pocket for the sample description	SK001.0010	1 pack-10 pcs
		SK001.0250	1 pack-250 pcs (25 x 10pcs)

MECCONTI S.A.R.L. SP. Z O.O.

T: +48 22 3974533, F: +48 22 5047213
sales@mecconti.com, support@mecconti.com
www.mecconti.com,
www.facebook.com/Mecconti
www.linkedin.com/company/mecconti