

Hy ConBOOTS



COMPLETE SETS DESIGNED FOR SAMPLING IN POULTRY HOUSES
TOWARDS SALMONELLA SPP. CONTAMINATION

SIMPLE TO USE - ALLOWS MEETING THE REQUIREMENTS IN EASIEST AND SAFEST WAY


COMPLETE SOLUTION - FROM THE MOMENT OF ENTRY INTO THE POULTRY HOUSE UNTIL
THE DELIVERY OF THE SAMPLE TO THE LABORATORY

COMPLIES WITH THE SAMPLING METHOD AS DESCRIBED IN EU COMMISSION
REGULATIONS (EU) NO 200/2010, (EU) NO 517/2011, (EU) NO 1190/2012


PRINCIPLES FOR SAMPLE COLLECTION

Faecal samples can be collected using sterile overshoes soaked in liquid e.g. MRD - Maximum Recovery Diluent. Depending on the type of flock, use the right number of sampling overshoes in one collection (for laying hens and turkeys - 2 pairs of sampling overshoes and for broiler flocks and breeding flocks - 5 pairs of sampling overshoes).


After putting on sterile sampling overshoes, you should go through the poultry house along the appropriate route, providing a representative sample of faeces. The collected samples should be gently slipped from the footwear and then placed in a sterile pouch, taking care that the collected faeces remain on the cladding. When putting on and taking off the sampling overshoes it is worth to use sterile gloves to prevent cross-contamination and transmission of contaminants from the sampling person's hands.

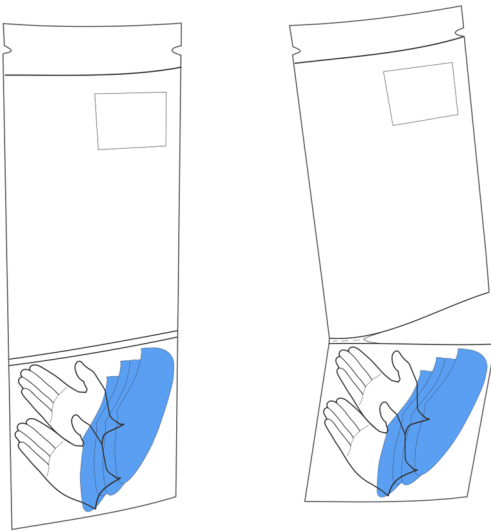


1. tear to separate

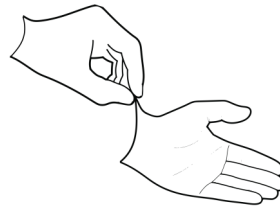


2. put the gloves on

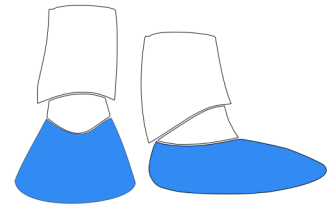

3. put sampling overshoes on

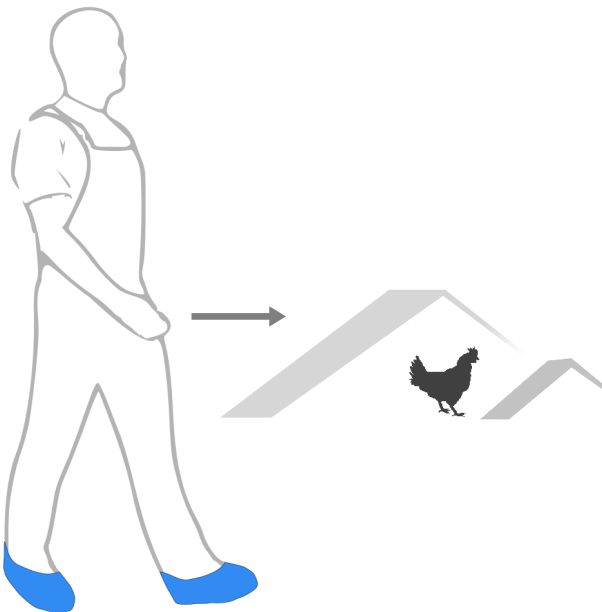

4. go through the poultry house along the appropriate route

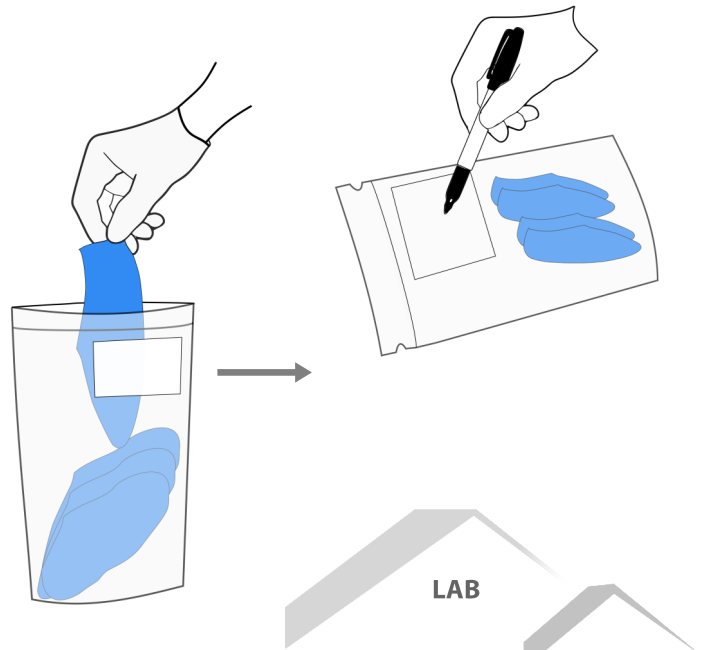

5. return used overshoes into sample bag



6. describe the sample and deliver to laboratory

HY CON BOOTS




- **2 OR 5 PAIRS SAMPLING OVERSHOES**
- **DISPOSABLE GLOVES IN ALL SETS**
- **EXTRA PROTECTION PLASTIC SHOE COVERS**
IN HY CON BOOTS PLUS SETS

SAMPLING OVERSHOES PRE-MOISTENED WITH MAXIMUM RECOVERY DILUENT

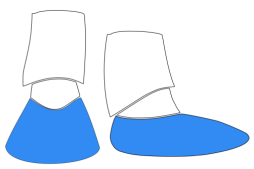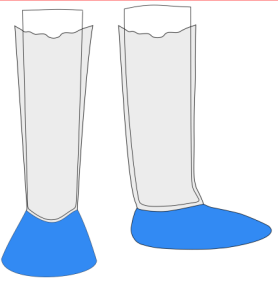		CODE	DESCRIPTION	ORDERING/PACKING	
	HY CON BOOTS	SK040	Hy Con Boots/MRD Sampling overshoes pre-moistened with MRD - 5 pairs Set include: pre-moistened sampling overshoes (5 pairs), 2 sample dual bags & 1 pair of disposable gloves	SK040.0025	1 pack - 25 pcs
		SK041	Hy Con Boots/MRD Sampling overshoes pre-moistened with MRD - 2 pairs Set include: pre-moistened sampling overshoes (2 pairs), sample dual bag & 1 pair of disposable gloves	SK041.0050	1 pack - 50 pcs
	HY CON BOOTS PLUS	SK042	Hy Con BootsPlus/MRD Sampling overshoes pre-moistened with MRD - 5 pairs Set include: pre-moistened sampling overshoes (5 pairs), 1 pair of extra protection plastic shoe covers, 2 sample dual bags & 1 pair of disposable gloves	SK042.0020	1 pack - 20 pcs
		SK043	Hy Con BootsPlus/MRD Sampling overshoes pre-moistened with MRD - 2 pairs Set include: pre-moistened sampling overshoes (2 pairs), 1 pair of extra protection plastic shoe covers, 1 sample dual bag & 1 pair of disposable gloves	SK043.0040	1 pack - 40 pcs

HyCon BOOTS MAP OF PRODUCTS

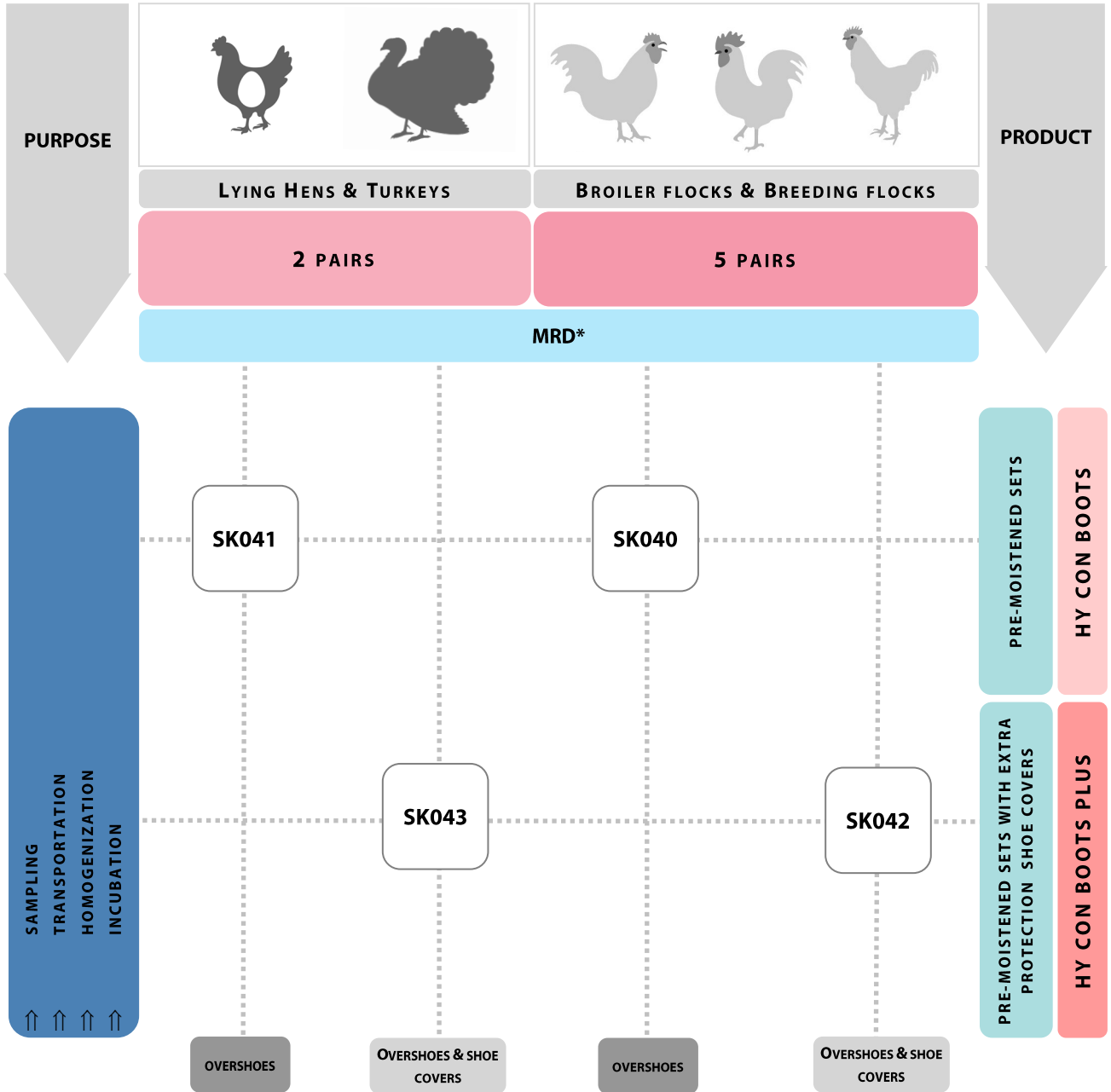

LEGENDS
 *MRD- Maximum Recovery Diluent

MECCONTI S.A.R.L. SP. Z O.O.

UL. RAKOWIECKA 36
02-532 WARSAW, POLAND

T: +48 22 3974533

F: +48 22 5047213

www.mecconti.com
sales@mecconti.com
www.facebook.com/Mecconti